

Microlink 751

User Manual

Biodata Limited

Manual Code: 751PDF-1.2

Issue Date: November 2015

Information in this document is subject to change without notice.
For updates see <http://www.microlink.co.uk/techsupp.html>

© Biodata Limited 2015
10 Stocks Street
Manchester M8 8QG
UK
Telephone: +44 (0)161 834 6688
Facsimile: +44 (0)161 833 2190
E-mail: sales@microlink.co.uk
<http://www.microlink.co.uk/>

Table of Contents

| | | |
|----------|---|------------|
| 1 | Introduction | 1.1 |
| 2 | Installing the Microlink 751 USB Unit | 2.1 |
| | 2.1 Introduction | 2.1 |
| | 2.2 Plugging the Microlink 751 into the Computer | 2.1 |
| | 2.3 The Microlink's Lights | 2.2 |
| | 2.4 Using Several Microlink 751 Units | 2.2 |
| | 2.5 Analogue Input Connection Notes | 2.4 |
| | 2.6 Digital Input and Output Connection Notes | 2.9 |
| | 2.7 Counter Connection Notes | 2.11 |
| | 2.8 Digital Input/Output and Counter Pin Connections Table | 2.12 |
| | 2.9 Specifications | 2.14 |
| | 2.10 The 59x Screw Terminal Units | 2.16 |
| 3 | Installing and Configuring Windmill Software | 3.1 |
| | 3.1 Introduction | 3.1 |
| | 3.2 What is Windmill Software? | 3.1 |
| | 3.3 Installing Windmill Software | 3.1 |
| | 3.4 Configuring Your System | 3.3 |
| 4 | Using Windmill Software with the Microlink 751 | 4.1 |
| | 4.1 Introduction | 4.1 |
| | 4.2 Options in SetupIML | 4.2 |
| | 4.3 The Rest of the Windmill Software Suite | 4.9 |

Introduction

Thank you for purchasing the Microlink 75I Measurement and Control Package. This Manual tells you:

- About the Microlink 751 unit and how to connect your signals (Chapter 2)
- How to install the Windmill software (Chapter 3)
- How to use Windmill software with the 751 unit (Chapter 4)

To use the 751 system you need a PC with a USB port. It should be running Windows 98SE, Windows 2000 or later.

Installing the Microlink 751 USB Unit

2.1 Introduction

The Microlink 751 unit plugs into the USB (universal serial bus) socket of your PC. It offers 16 analogue inputs, 32 digital inputs/outputs and 8 counters.

The 751 has two 37-way connectors: 1 for analogue connections and one for digital and counter connections. If you prefer you can make your connections to screw terminals, using one of the optional Microlink 59x range of external units. You will also need a 59x unit if you plan to interface transducers like thermocouples or strain gauge bridges.

This manual tells you how to connect your signals to the Microlink 751, and about the 59x screw terminal units. After making your connections see the next chapter for details of installing the Windmill software.

2.2 Plugging the Microlink 751 into the Computer

Plugging the Microlink into your PC could not be easier: just use the USB cable provided. You don't have to switch off your computer first—or even restart Windows.



The Microlink 751 is powered from the USB port: you don't need an extra power supply box. If you have several 752 units, make sure their ID codes are set (Section 2.4) before plugging them into the PC.

2.3 The Microlink's Lights

The Microlink 751 has two red lights labelled ENUM and BUSY. Neither of these will come on until you have installed the Windmill software.

- ENUM ENUM stands for Enumerated. This is lit when the Microlink has been powered on by the USB plug and play controller. It is a good indication that the Windmill USB driver software has been correctly loaded.
- BUSY This is lit for the duration of each USB communication. It is not active until the Microlink has been enumerated.

2.4 Using Several Microlink 751 Units

As you can connect up to eight Microlink 75x units to the USB, each one must have some way of identifying itself. This is provided by 3 pins on the digital connector. (If you have just the one Microlink 751, its ID Code will be 0 and you can ignore this section.)

The pins in question are numbers 28, 9 and 10. You set an ID code by connecting these pins to 0 V, as follows.

| ID Code | Pin 10 | Pin 9 | Pin 28 |
|---------|--------|-------|--------|
| 0 | no | no | no |
| 1 | no | no | yes |
| 2 | no | yes | no |
| 3 | no | yes | yes |
| 4 | yes | no | no |
| 5 | yes | no | yes |
| 6 | yes | yes | no |
| 7 | yes | yes | yes |

See the Pin Connections Table on page 2.8.

Make a note of the ID Codes, you'll need them when using Windmill ConfIML to install the driver software. You may find it helpful to label your 751s with their codes—for quick identification.

Set the ID code before connecting the Microlinks to the PC.

2.5 Analogue Input Connection Notes

The Microlink 751 provides 16 differential analogue inputs. It also has an auxiliary channel which is reserved for special inputs such as cold junction measurement or bridge excitation monitoring (used with thermocouples and strain gauges). It uses an integrating analogue-to-digital converter, where the integration time and resolution are under software control (set in the Windmill ConfIML program, which starts automatically when you install the software).

At regular intervals Windmill uses a stable on-board reference voltage for recalibration. Use ConfIML to set this recalibration interval.

2.5.1 Differential Inputs

All the inputs are differential: for each input signal there are two signal wires. The measurement is the difference in voltage between the two wires. The two signals go into separate high-impedance amplifiers which monitor the voltage between the input and ground. The outputs of the two amplifiers are then subtracted to give the difference between the + and – inputs. For small signals differential inputs are much better than single-ended inputs, because the subtraction of the voltages on each of the input wires means that any voltage common to both wires is removed, so reducing noise.

It is important to remember that the voltages at both inputs must be within the amplifier operating range. A classic error is to connect a battery between positive and negative inputs with no other connection. Although the difference between the inputs is well defined, the actual voltage at each input could be anything. Connecting one end of the battery to the 0 V input, either direct or via a resistor, would solve the problem.

The 0 V of the Microlink 751 is connected to computer earth, which is usually connected to the mains earth. When making your connections you should follow this policy.

- a. If your signal source is “floating”, i.e. has no reference to mains earth, then you must provide a reference by connecting one end of

it to the 0 V input, either direct or via a resistor. The resistor could have any value up to several M Ω . However large values could cause 50 Hz problems if your signal source has much leakage to earth.

- b. If your signal source is itself earthed then you should connect only positive and negative inputs. You should make no connection to Microlink 751 earth.

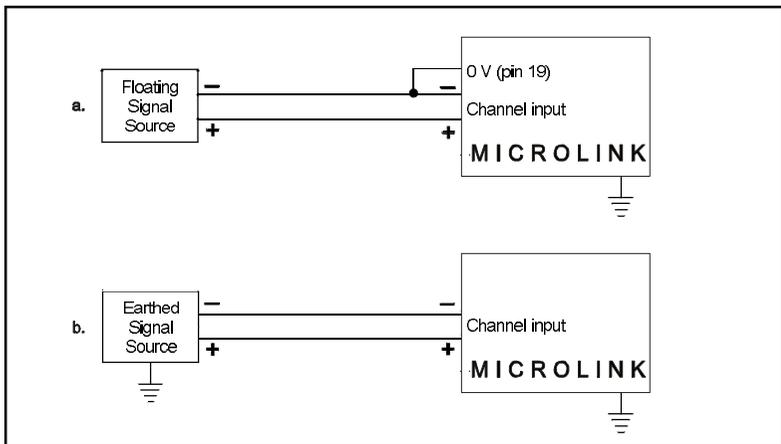


Figure 2.1 For floating signals you need to connect one end of the signal to the 0 V input on pin 19

2.5.2 Input Voltage Range

The Microlink 751 operates correctly with input voltages in the range ± 11 V. The inputs will reject voltages which are common to both positive and negative inputs. These common mode voltages could be as big as 13 V.

2.5.3 Maximum Input Voltage

The input multiplexers are protected against dc voltages of 33 V above the power supply. This means ± 48 V if the computer is switched

on, ± 33 V if it is switched off. If the over-voltage is transient then protection extends as far as ± 300 V. When a voltage above the power supply is applied to the unit its protection mechanism comes into action, and this draws some current from the signal source. This effect can be a problem when the computer is switched off as it now draws current from any signal. This current is limited by 4K7 resistors. Extra series resistors can be added to reduce this fault current.

2.5.4 Unconnected Inputs

You can leave unused inputs unconnected, but if you attempt to read from these unconnected inputs do not expect to get 0 V. They could be any value. If another connected channel has recently been read, the unconnected input will return a similar value. This is not crosstalk. It occurs because the input capacitance of the amplifier is charged to the voltage of the previous channel and has little incentive to change when connected to an open circuit.

2.5.5 Auxiliary Input

This input has all the facilities of the other sixteen. In Windmill software it is reserved for measuring cold junction temperatures with thermocouples or excitation voltage in bridge circuits. If using a 593 thermocouple input unit or a 594 bridge input unit, the auxiliary input is automatically connected as required by the software.

2.5.6 Use with 593 Thermocouple Input Unit

When used with this unit the auxiliary input is automatically connected to the cold junction sensor. This is an RTD (resistance temperature device) with 0.8 milliampere flowing through it. If inputs are unconnected the thermocouple break detection circuitry will cause them to read positive full scale.

Pins 18 and 37 (-15 V and $+15$ V) provide a power supply for the 593—don't connect any other signals to these pins.

2.5.7 Use with 594 Bridge Input Unit

When used with this unit the auxiliary input automatically measures the excitation voltage connected to the 594.

2.5.8 Analogue Input Pin Numbers

Make your connections to the analogue 37-way D connector as detailed in the table on page 2.8.

Microlink 751 - 16 Analogue Inputs

| | | | |
|-------------|----|----|-------------|
| +15 V | 37 | 19 | 0 V |
| + Auxiliary | 36 | 18 | -15 V |
| + Input 15 | 35 | 17 | - Auxiliary |
| + Input 14 | 34 | 16 | - Input 15 |
| + Input 13 | 33 | 15 | - Input 14 |
| + Input 12 | 30 | 14 | - Input 13 |
| + Input 11 | 31 | 13 | - Input 12 |
| + Input 10 | 30 | 12 | - Input 11 |
| + Input 9 | 29 | 11 | - Input 10 |
| + Input 8 | 28 | 10 | - Input 9 |
| + Input 7 | 27 | 9 | - Input 8 |
| + Input 6 | 26 | 8 | - Input 7 |
| + Input 5 | 25 | 7 | - Input 6 |
| + Input 4 | 24 | 6 | - Input 5 |
| + Input 3 | 23 | 5 | - Input 4 |
| + Input 2 | 22 | 4 | - Input 3 |
| + Input 1 | 21 | 3 | - Input 2 |
| + Input 0 | 20 | 2 | - Input 1 |
| | | 1 | - Input 0 |

Please read the Connection Notes on the previous pages before making your connections.

2.6 Digital Input and Output Connection Notes

The Microlink 751 provides digital input to the computer and output control by the computer. Its 32 general purpose input and output lines are arranged in 4 groups or ports. Each port can be either input or output (set using the Windmill SetupIML program). All ports power-up as inputs. The ports are referred to as Port 0 to Port 3. Port 3 also functions as 8 event counters, detailed in the next section.

2.6.1 Input Voltages

All inputs are high impedance CMOS type. They are TTL and 5 V CMOS compatible. Input Voltages should be within the range 0 to 5 V. Higher Voltages can be dealt with by the addition of resistor networks. This can be conveniently done on a 590 unit. Input protection can be provided in a similar manner.

2.6.2 Contact Closures

You can interface to contact closures using a resistor to tie the input to either 5 or 0 V. The contact then switches the line to either 0 or 5 V. The resistor can be fitted to a 590 unit.

2.6.3 Noisy Inputs

Input Filters can be fitted to a 590 unit if required.

2.6.4 Output Drive

The outputs are TTL and 5 V CMOS compatible. They can drive 15 LSTTL loads. You can increase the output drive by using additional transistors, which can be fitted to the 590 unit. Currents of 1 amp can easily be switched.

2.6.5 Power-Up State

The 751 unit will power-up as all inputs. If you intend to use the card to control outputs then you may want to define logic states at power-up. This can be done by resistors which tie the lines to either 0 or 5 V, mounted on a 590 unit.

2.6.6 Pin Numbers

Make the I/O connections to the digital 37-way connector. See the Pin Connections Table on page 2.13.

2.7 Counter Connection Notes

The 751 provides eight 16-bit totalise (event) counters which can each count up to 65535. These are located on Port 3 of the digital I/O connector. If you are using counters, set Port 3 as an input only, using the Windmill SetupIML software.

The 751 unit monitors the state of the 8 input lines once every millisecond and maintains a count for each of them. It does this whether or not you intend to use the lines as counters. You can still read Port 3 as a normal digital input, even if you are also using it to count.

2.7.1 Input Voltages

See Section 6.12.6.1 for safe voltage levels.

2.7.2 Count Inputs

A valid count is declared if the input is low for 2 milliseconds then high for two milliseconds. This gives a theoretical maximum count speed of 250 Hz.

2.7.3 Pin Numbers

Make the counter connections to Port 3 of the Digital Connector. The counter pin numbers are on page 2.13.

2.8 Digital Input/Output and Counter Pin Connections Table

The pin numbers given on the next page are those for the digital 37-way D socket.

The analogue input pin numbers are on page 2.8.

*Microlink 751 - Pin Connections for
Digital Inputs/Outputs and Counters*

| | | | |
|--------------|----|----|--------------|
| | | 19 | 0V |
| Port 3 Bit 0 | 37 | 18 | Port 3 Bit 1 |
| Port 3 Bit 2 | 36 | 17 | Port 3 Bit 3 |
| Port 3 Bit 4 | 35 | 16 | Port 3 Bit 5 |
| Port 3 Bit 6 | 34 | 15 | Port 3 Bit 7 |
| Port 2 Bit 0 | 33 | 14 | Port 2 Bit 1 |
| Port 2 Bit 2 | 32 | 13 | Port 2 Bit 3 |
| Port 2 Bit 4 | 31 | 12 | Port 2 Bit 5 |
| Port 2 Bit 6 | 30 | 11 | Port 2 Bit 7 |
| Not used | 29 | 10 | ID Code 2 |
| ID Code 0 | 28 | 9 | ID Code 1 |
| Port 1 Bit 0 | 27 | 8 | Port 1 Bit 1 |
| Port 1 Bit 2 | 26 | 7 | Port 1 Bit 3 |
| Port 1 Bit 4 | 25 | 6 | Port 1 Bit 5 |
| Port 1 Bit 6 | 24 | 5 | Port 1 Bit 7 |
| Port 0 Bit 0 | 23 | 4 | Port 0 Bit 1 |
| Port 0 Bit 2 | 22 | 3 | Port 0 Bit 3 |
| Port 0 Bit 4 | 21 | 2 | Port 0 Bit 5 |
| Port 0 Bit 6 | 20 | 1 | Port 0 Bit 7 |

•

Please read the Connection Notes on the previous pages before making your connections.

2.9 Specifications

2.9.1 Microlink 751 Unit

| | |
|--|----------------|
| Dimensions (mm) | 180 x 120 x 40 |
| Maximum number of 751s | 8 |
| Maximum length of cable | 5 m per cable |
| Maximum distance from PC can be increased by use of USB hubs | |

2.9.2 Analogue Inputs

| | |
|-------------------------------------|--------------------------------------|
| Number of inputs | 16 |
| Maximum safe input voltage | |
| Computer on | ±48 V |
| Computer off | ±33 V |
| Transient | ±300 V |
| Ranges (set from software) | ±10 V, ±1 V, ±100 mV, ±10 mV |
| Common mode range | ±13 V |
| Relative accuracy of ranges | |
| gain = 1000 | ±0.1% |
| gain = 1, 10, 100 | ±0.05% |
| A-D performance (set from software) | |
| Resolution (bits) | Integration Time (msec) |
| 12 | 2.5 |
| 13 | 5 |
| 14 | 10 |
| 15 | 20 |
| 16 | 40 |
| 18 | 160 |
| Maximum speed of sampling | 10 samples per second per channel |
| up to a maximum of | 80 samples per second |
| Maximum linearity error | ±0.02% |
| Input impedance | 100 MΩ |

2.9.3 Digital Inputs and Outputs

| | |
|---------------------------|-------------------------|
| Maximum number of inputs | 32 |
| Maximum number of outputs | 32 |
| Power-up state | all inputs |
| Maximum speed | 160 channels per second |
| Voltage Inputs | |
| Compatibility | TTL and 5 V CMOS |
| Range | 0 to 5 V |
| Voltage Outputs | |
| Compatibility | TTL and 5 V CMOS |
| Drive | 15 LSTTL loads |

2.9.4 Counters

| | |
|----------------------------|-----------------------|
| Maximum number of counters | 8 |
| Resolution | 16 bits |
| Type of counters | Event (totalise) |
| Maximum count speed | 160 counts per second |
| Compatibility | TTL and 5 V CMOS |
| Input voltage range | 0 to 5 V |

2.10 The 59x Screw Terminal Units

The Microlink 59x range provides screw terminals for the Microlink 751, allowing signals to be permanently wired. These boxed 59x units sit next to the Microlink 751. Some of the 59x range allow transducers like thermocouples and strain gauge bridges to be monitored by the 751. Extra facilities are available when components are fitted, such as protection from high voltages. There are 4 units in the range.

| | |
|--------|---|
| 590 | Screw Terminals |
| 590-8A | Screw Terminals and Current Measurement |
| 593 | Screw Terminals and Temperature Measurement |
| 594 | Screw Terminals and Strain Measurement |

The 59x range is optional, so you can ignore the rest of this chapter if you have bought just the Microlink 751 and Windmill package.

2.10.1 590 Unit—Screw Terminals

The 590 is a boxed screw terminal card. Extra facilities are available with this card, including:

- noise filtering,
- input protection from high voltages,
- higher voltage inputs than normal,
- contact closure inputs,
- ac inputs,
- biasing to ground,
- external current source switching,
- programmable resistance,
- special output ranges.

Connection

The 590 unit has space for 3 rows of 18 screw terminals. The row nearest the front panel is connected to pin 19 of the analogue D connector on the front of the Microlink. This pin is connected to 0 V. The other 2 rows are connected to pins 1–18 and 20–37 as indicated by the small figures by the terminals.

Fitting Application Components

Series Components

Component positions S1 to S37 (0.4 inch pitch) are available to mount components in series with D-type pins 1 to 37. Only S1 and S37 are named on the PCB because of lack of space. However the numbering sequence is the same as that of the nearby B1 to B37 components. To fit a component in series with, for example, pin 10

- Place the component in the S10 position (level with B10),
- Solder it on the copper side,
- Cut the track under S10 at the point where it narrows.

The main application of series components is series protection resistors. These could have any value from hundreds of ohms to megohms depending on requirements.

Parallel Components

Component positions P0 to P17 (0.2, 0.3, 0.4 inch pitch) are provided to mount components across analogue inputs 0 to 15, 16 is the auxiliary input, 17 completes the pattern. Components can be simply soldered into these positions.

Input Filters

Low pass analogue input filters can be produced by fitting resistors in the 2 series elements of an analogue channel, and a capacitor in the parallel element. For instance a filter can be implemented on channel 1 by fitting resistors to S2 and S21 and a capacitor to P1. Typical values would be 10 K Ω and 1 μ F. These would produce a cut off frequency of about 8 Hz.

2.10.2 590-8a Unit—Screw Terminals with Current Measurement

The 590-8A is a boxed milliamp current detection card, arranged as 8 current and 8 voltage inputs. The same extra facilities are available with this card as with the 590.

2.10.3 593 Unit—Screw Terminals with Temperature Measurement

The 593 is an isothermal box specialised for thermocouple monitoring of temperatures. It includes a cold junction sensing RTD with an associated constant current source (0.8 milliamps). The RTD determines the temperature at the termination point. This reference temperature is used with the analogue input channel's voltage reading to calculate the temperature at the thermocouple junction.

The Windmill software automatically connects the cold junction to the auxiliary channel and compensates for it. It reads the channel directly, showing the cold junction temperature in Celsius.

Protection from high voltage inputs is provided when series resistors are added. Noise problems can be solved by fitting an input filter and broken thermocouple leads can be detected.

The cold junction sensing circuitry is located between the two leftmost rows of screw terminals. This circuitry is connected by the PCB to the auxiliary channel. Bias resistors are fitted to produce a small current through the thermocouples. This allows thermocouple break detection.

2.10.4 594 Unit—Screw Terminals with Strain Measurement

The 594 is a boxed 16 bridge inputs card, which enables the 751 unit to monitor strain gauge bridges and balanced bridges such as pressure transducers.

For strain gauges you need an external excitation voltage, which can supply sufficient current to keep all the bridge circuits energised. Four sets of 16 screw terminals give + and – excitation and + and – signals for each of 16 bridges. Two 1 k Ω termination resistors are mounted in half bridge configuration whilst high quality 350 and 120 Ω resistors can be provided for the completion of quarter bridges. The auxiliary channel of the 751 unit monitors the excitation voltage. The A-D converter is suitable for direct measurement of voltage imbalance provided you choose a high resolution.

When monitoring a bridge input, the Windmill software automatically reads excitation voltage and performs the bridge calculation to produce a reading in microstrain. You can set a zero reference level and monitor changes relative to that level.


With the 594 unit you can configure each input channel for any of the following inputs.

| | |
|----------------|---|
| voltage input | |
| quarter bridge | single strain gauge |
| half bridge | tensile + compressive strain gauge |
| half bridge | normal + transverse strain gauge |
| full bridge | 2 tensile + 2 compressive gauges |
| full bridge | 2 normal + 2 transverse gauges |
| full bridge | tensile normal + compressive normal + tensile transverse + compressive transverse gauges |

The figure on page shows the general arrangement of the 594. For each input channel (0–15) there are 5 screw terminals. These are arranged in 5 rows labelled, 0V, +EX, –EX, +IN, –IN, with channel number printed by each terminal. Common to all channels are the

POWER INPUT terminals and two precision resistors connected across the power supply to form a half bridge. The power input terminals are directly connected to the auxiliary channel of the 751. Windmill software reads this channel to measure the excitation voltage for use in the bridge equations. Each channel has associated with it 2 link mounting positions and one resistor position. These are:


- RX Position for mounting a quarter bridge completion resistor
- QX Link to short out the quarter bridge completion resistor
- HX Link to connect the half bridge to the -IN of channel X

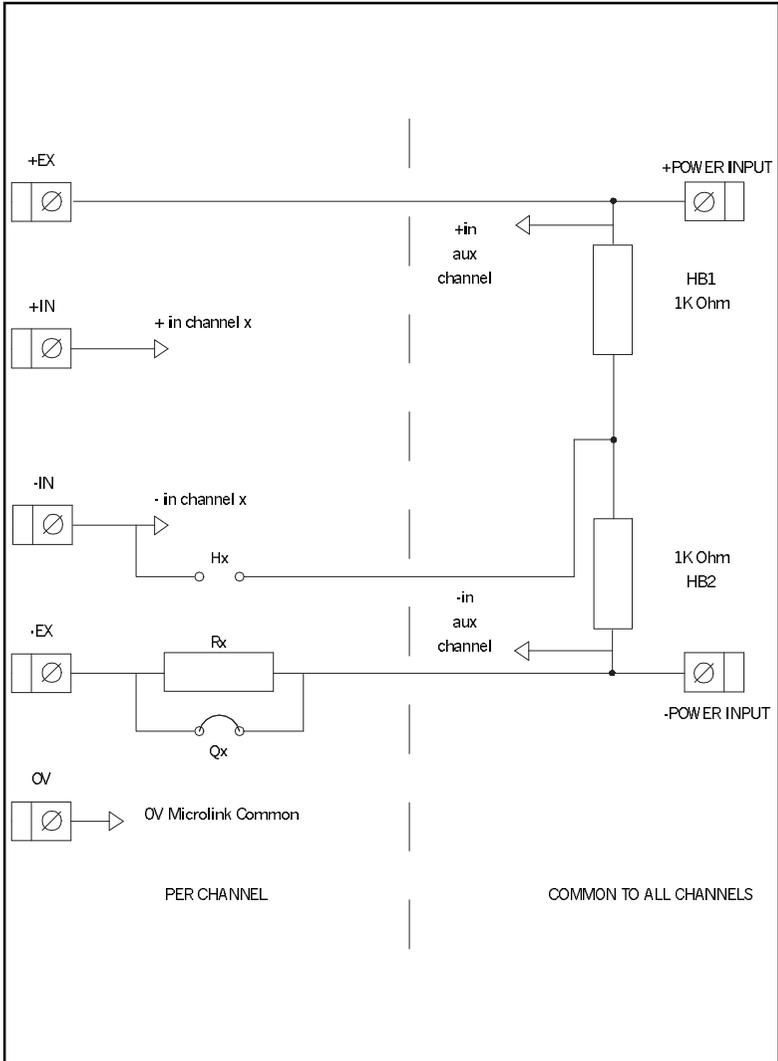

General Arrangement of 594

Full Bridge Connection


The figure below shows a full bridge connection. The QX quarter bridge link must be fitted to connect the –POWER INPUT to the –EX. Four wires are needed to connect your bridge and the earth is available for a cable screen.

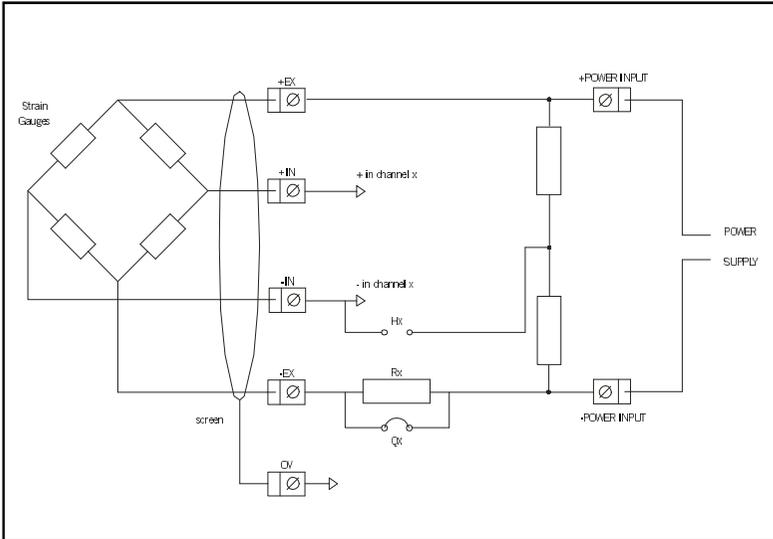

594 Full Bridge Connection

Half Bridge Connection

This figure shows a half bridge connection. The HX and QX links must both be fitted. The HX link connects the half bridge to the –IN of the channel, whilst the QX link connects the –POWER INPUT to the –EX. Three wires are needed to connect your bridge and the earth is available for a cable screen.

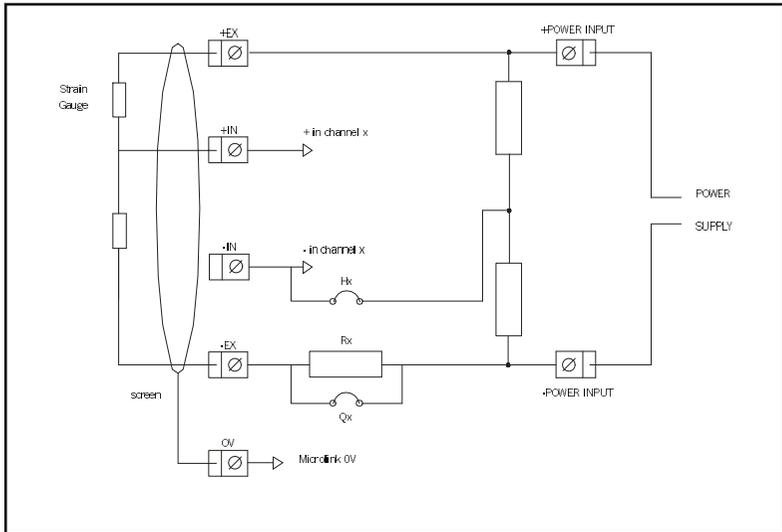


594 Half Bridge Connection

Quarter Bridge Connection

The figure below shows a quarter bridge connection. The HX link connects the half bridge to the -IN. The RX bridge completion resistor must be fitted. This resistor should be a high precision, low temperature coefficient device. Its value should be the same as the nominal resistance of your strain gauge, 120 and 350 Ω being the commonest values. In other applications it should be chosen to balance the bridge near the centre of your measurement range. For instance if you are making precise RTD measurements around ambient temperature then 100 Ω would be suitable. If however you were monitoring higher temperatures a suitably larger value would be chosen.

It is normal practice when making a quarter bridge connection to take all 3 wires to the transducer. This balances out the effects of lead resistance.

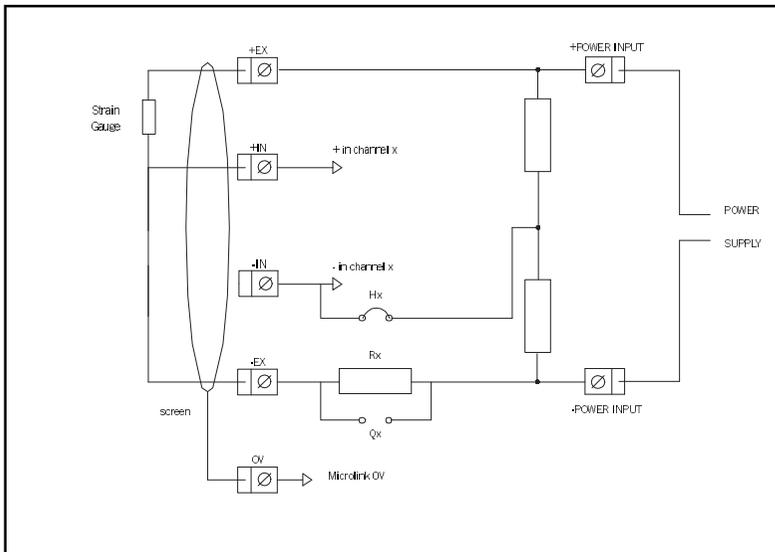

594 Quarter Bridge Connection

Installing and Configuring Windmill Software

3.1 Introduction

There are two steps to installing and configuring the Windmill measurement and control software.

1. Install the software onto the hard disk.
2. Tell the software about your measurement hardware.

For installation you will need a personal computer with a USB port that is running Windows 98SE or later.

3.2 What is Windmill Software?

Windmill is a ready-to-run suite of applications for data acquisition and control. You can be up and running in very little time as no programming is required. The standard suite includes data logging, charting and output control applications. You can also transfer data directly into third-party applications like Microsoft Excel or Access. Other Windmill modules are available—see the Windmill Software Internet catalogue for details <http://www.windmillsoft.com/>.

3.3 Installing Windmill Software

1. Connect the 751 data acquisition unit to your computer, as detailed in Chapter 2.
2. Install the Windmill software:
Insert the CD. The Installation software should run automatically. If not select Run from the Start menu and type d:setup (where d: is your CD drive).

3. When installation is complete the Windmill Configuration program, ConfIML, automatically starts. This asks for details of your hardware. See Section 3.4 for details.

3.3.1 Upgrading From Earlier Versions of Windmill

You can run Windmill alongside earlier versions of the software. Once you are happy that you no longer need a previous copy of Windmill simply run its “Uninstall” program to delete it.

3.3.2 Running Several Copies of Windmill

Unless you have bought a multiple licence, you are only permitted to run one copy of the software at any one time. Please contact your supplier if you require a multiple licence.

3.3.3 Installing under Windows 7, 8 and Later

If you are using Windows 8 or 10, and you encounter problems with your USB devices, see <http://www.windmill.co.uk/usb-device-windows8.html>

3.4 Configuring Your System

The Windmill Configuration program, ConfIML, records the details of your data acquisition hardware. It will run at the end of the installation process, and you should run it again from Windows whenever your acquisition hardware changes—for example when you install additional units.

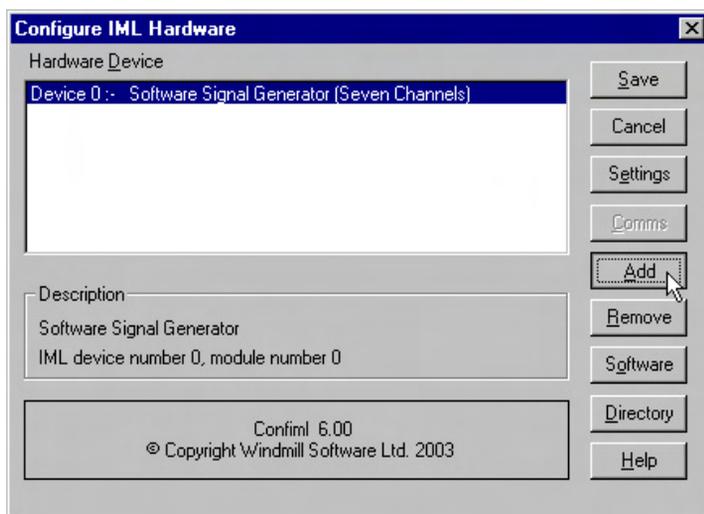


Figure 3.1 The ConfIML Main Window

The first thing you need to do is press the Add button to include your Microlink 751 unit.

3.4.1 Adding New Hardware

The Add IML Hardware dialogue lists the acquisition and control devices for which you have installed drivers. Select the Microlink 751 unit and press the Add button. This will take you to the Hardware Settings dialogue (Section 3.4.2).

3.4.2 Changing Hardware Settings

ConfIML needs to know some information about your 751 unit. Press the Help button or see below if in doubt about the answers.

Which Type of Card do you have?

Choose the Microlink 751.

ID Code of the 751 Unit?

As you can connect several 751 units to your PC, you need to tell ConfIML which one you are currently configuring. Do this by selecting the correct ID Code. If you have only one 751 its ID Code is 0. (There should be as many ID codes as there are 751 units connected, and each code should be different. If not, change the codes on individual units as detailed in Section 2.2.4.)

Which Type of Transducer Connection Unit is Attached?

There are three options. Which you choose depends on the 59x units you have connected.

None/General Purpose

Choose this if you:

- are not using a 59x unit,
- are using a 590 screw terminal voltage unit, or
- are using a 590-8A screw terminal current unit.

Thermocouple

Choose this if you are using a 593 thermocouple terminal unit.

This monitors the temperature of connections, which is essential when using thermocouples. You can use channels not required for thermocouples for general voltage measurement.

Strain Bridges

Choose this if you are using a 594 strain gauge bridge connection unit.

This monitors an external bridge excitation supply so that

strain gauge outputs can be converted into readings in microstrain. Again, you can use channels not required for strain for general voltage measurement.

Select the Integration Time (Resolution)

This allows you to select the trade off between speed, precision and noise rejection. Seven settings are available—the slower the speed, the better the precision and the more random noise is averaged out. The exact speeds and resolutions vary with the computer, transducer type and so on, but the following table gives some typical figures.

| Setting | Typical Speed | Voltage Resolution |
|---------|---------------|----------------------|
| 12-bit | 80 reads/sec | 5 mV in ± 10 V |
| 13-bit | 64 reads/sec | 3 mV in ± 10 V |
| 14-bit | 48 reads/sec | 1.5 mV in ± 10 V |
| 15-bit | 32 reads/sec | 0.8 mV in ± 10 V |
| 16-bit | 16 reads/sec | 0.4 mV in ± 10 V |
| 18-bit | 6 reads/sec | 0.1 mV in ± 10 V |

Settings from 15-bit to 18-bit will be effective at removing 50 Hz mains noise as these integrate over one or more complete mains cycles.

The resolution figures show the ability to detect small changes. The absolute accuracy will depend on several factors including your transducers.

Interval Between Automatic Recalibrations (in Minutes)?

This selects how frequently the software pauses to perform a recalibration of the Microlink 751 using its stable on-board reference voltages. This will cancel any errors introduced by temperature or power supply changes since the last recalibration, but can take up to a second to complete. During this time, no readings may be taken. You can vary the interval from 1 to 1440 minutes (24 hours), but every 20 minutes is generally sufficient.

3.4.3 751 Devices

After making your choices and pressing OK, the ConfIML Summary window appears showing your 751 as three hardware devices: 751 analogue inputs, 751 digital ports and 751 event counters.

3.4.4 Setting the Working Directory

When you use Windmill you will create two types of files, those which hold data and those which hold the Windmill programs' settings. You can specify a default folder in which to store these files— known as the Working Directory. Initially the working directory is set to wherever you installed Windmill. To change this, press the Directory button.

Choose a new directory (folder) and press Save. Windmill will create two sub-directories under this choice, called setup and data. By default, all the Windmill settings will be stored in “setup” and all the data you collect stored in “data”.

3.4.5 Saving the Settings

Save your settings and they will be used every time you run Windmill. You do not need to run ConfIML again, unless you add more units or want to change the settings.

After closing ConfIML start the SetupIML program and choose how you wish to use the hardware. Refer to the next chapter and SetupIML's Help file (setupiml.hlp) for details.

3.4.6 The Software Signal Generator

In addition to data acquisition and control hardware, ConfIML lists the Software Signal Generator. This is a special driver which simulates a device with seven channels, each channel producing a different signal. No special hardware is required—the data values are produced by calculation.

The Software Signal Generator lets you experiment and practise with Windmill, without being concerned about the hardware

To install the signal generator you first need to add it to the list of devices. In the ConflIML Summary window press the Add button (Section 3.4.1). Select Software Signal Generator and again press the Add button. You're taken to the Hardware Settings dialogue (Section 3.4.2) where you can choose options for five of the channels. Press the Help button in this dialogue for more information on the signal generator.

3.4.7 The IML Device Icon

Whenever you run a Windmill program one or more IML Device icons will appear. Different icons identify different hardware drivers, software signal generators and so on. The Windmill applications can't run without these, so don't close them whilst using Windmill.

Using Windmill Software with the Microlink 751

4.1 Introduction

This chapter explains which settings in the Windmill SetupIML program apply to the 751, and about the rest of the Windmill software suite.

You can use the Microlink 751 to monitor voltages. With additional Microlink 59x units, you can also monitor currents, thermocouples and strain gauge bridges. If you wish to measure resistance or pH, please contact Windmill Software for advice.

Before starting Windmill make sure your 751 is plugged into the computer's USB port, and, if you are using a 59x unit, that it is plugged into the 751's analogue connector.

4.2 Options in SetupIML

SetupIML is the Windmill program that lets you save libraries of setup files, each holding details about how you want to use individual channels. A Windmill 751 is shown as three devices in SetupIML: Analogue Inputs, Digital Inputs and Event Counters. Choose one from SetupIML's Device menu. Now double-click a channel to configure it. Full details of using SetupIML are in its Help file (setupiml.hlp) and Chapter 3 of the *Windmill User Manual*.

4.2.1 General Purpose Analogue Input Connector

The input channels are numbered 0 to 15. Using SetupIML you can configure each channel as follows:

- enable or disable
- re-name
- set to a specific input range or allowing the software to pick the range automatically
- give a new units name, scale factor and offset
- set alarm levels

4.2.2 Thermocouple Input Connector

The signal inputs are numbered 0 to 15. Using SetupIML you can configure each channel as follows:

- enable or disable
- choose one of six common thermocouple types for automatic linearisation to Celsius, or choosing to read raw voltage
- re-name
- set to a fixed range or allowing the software to pick the range automatically
- give a new units name, scale factor and offset
- set alarm levels.

Channel 16 is normally disabled, but if enabled you can use it to read the cold junction reference temperature at the screw terminals. This can be used to linearise the voltages from other types of thermocouple in some external software such as a spreadsheet.

4.2.3 Strain Gauge Bridge Connector

The signal inputs are numbered 0 to 15. Using SetupIML you can configure each channel as follows:

- enable or disable
- set to one of six bridge configurations (see below) or set to read raw voltage
- re-name
- set to a fixed range or allowing the software to pick the range automatically
- set to a gauge factor
- if transverse strain is being measured, specify Poisson's ratio for the material under test
- set alarm levels.

See also the Special commands described on page 4.5.

The supported bridge configurations are:

Quarter bridge

One active gauge and three fixed resistors.

Half bridge

One gauge measuring tensile strain (+E), one measuring compressive strain (-E), and two fixed resistors.

Half bridge

One gauge measuring normal strain (+E), one measuring transverse strain (-vE), and two fixed resistors. Poisson's ratio must be specified.

Full bridge

Two gauges measuring tensile strain (+E) and two measuring compressive strain (-E).

Full bridge

Two gauges measuring normal strain (+E) and two measuring transverse strain (-vE). Poisson's ratio must be specified.

Full bridge

One gauge measuring normal tensile strain (+E), one measuring normal compressive strain (-E), one transverse tensile strain (+vE), and one compressive transverse strain (-vE). Poisson's ratio must be specified.

In all these cases the answer is reported directly in microstrain.

Channel 16 is normally disabled but if enabled it can be used to read the bridge excitation voltage, which must be the same for all channels, and is used in the calculations.

4.2.4 Special Hardware Commands

The channel dialogue boxes will have the “Special” button enabled. This allows access to the following commands:

Recalibrate

Forces an immediate recalibration of the whole board. This is in addition to the periodic calibration cycles which happen automatically.

Cal_Interval

Displays the time between automatic calibration cycles, as selected in ConflML (described in Section .

Resolution

Displays the resolution and integration time for this card, as selected in ConflML (described Section).

Show_Cal_Data

Used only during manufacturer's test procedures.

When a strain gauge bridge is in use, three additional commands are available:

Set_Zero_Now

This reads the bridge output ratio and uses it as the zero strain reference level. All subsequent readings will be changes in strain relative to that point, until another special command is issued or the software is closed down.

Describe_Offset

This shows the bridge output level which is currently being used as the zero strain reference level (as a fraction of excitation voltage).

Disable_Offset

Resets the zero strain reference to zero volts, so absolute readings are obtained. This is the default state each time the software starts running.

4.2.5 Digital Ports

The Microlink 751 unit provides four digital ports, each with 8 lines. You can use each port for input or output. You can explicitly control the choice, or it can be deduced automatically by the software. All ports start as inputs, but if any data is sent to a port by a Windmill program it immediately switches to output mode. Output lines can also be controlled by the alarm detection features of the input channels. This will cause the whole port to switch to output mode.

You can also use port 3 for event counting—in which case you must set it as an input.

The channels are numbered as follows:

| | |
|------|-------------------------------------|
| 0100 | Line 0 of port 0 |
| 0101 | Line 1 of port 0 |
| 0102 | Line 2 of port 0 |
| .. | ... |
| 0107 | Line 7 of port 0 |
| 0108 | Direction control signal for port 0 |
| 0109 | not used |
| 0110 | Line 0 of port 1 |
| 0111 | Line 1 of port 1 |
| ... | ... |
| 0117 | Line 7 of port 1 |
| 0118 | Direction control signal for port 1 |
| 0119 | not used |
| 0120 | Line 0 of port 2 |
| etc | |

In SetupIML you can replace the numbers with **meaningful names**.

By default, each channel is linked to a single line and the two digital states are called ON and OFF. You can change these names to more suitable ones, for example FAST and SLOW or OPEN and SHUT.

SetupIML lets you group the lines within each port into multi-bit channels which are displayed or controlled as single values transferred

via the first channel in the group. Multi-bit values can be chosen as binary, decimal or hexadecimal, e.g.

| Binary | Decimal | Hexadecimal |
|----------|---------|-------------|
| 11 | 3 | 3 |
| 1010 | 10 | A |
| 10001 | 17 | 11 |
| 11111111 | 255 | FF |

Channels 0108, 0118, 0128 and 0138 are the direction control signals for ports 0 to 3. When set to 0, the whole port is used for input; when 1, the port is used for output. These channels can be viewed and also altered, but they are normally disabled and must be explicitly enabled from SetupIML before use. As noted above, direction control can usually be sensed automatically by the software.

Channel 0109, 0119 and 0129 are not used and cannot be enabled.

4.2.6 Event Counters

The Microlink 751 unit provides eight 16-bit event counters. The default names are 0200, 0201 to 0207.

Each counter starts at zero and counts pulses on the corresponding input lines, to a maximum of 65535. They can be reset to zero at any time by sending "0" to the channel using the AnalogOut, Graphics or Test-Seq programs. No other values can be sent to the counters.

Use SetupIML to change the names, enable or disable each channel, and choose one of two operating modes—accumulating count or resetting count.

Accumulating Count

Simply keeps counting until explicitly reset.

Resetting Count

Starts again from zero after each reading. This shows the number of pulses since the last reading, but it can only be used where a single program is reading the counter. If several different programs were accessing the counter simultaneously, they would all be resetting it at different times, so the results would be unpredictable.

SetupIML also lets you apply a scale factor and offset to the count. For example, if the pulses came from a flow meter which produced one pulse for each 50 millilitres, then a scale factor of 0.05 would give a reading in litres.

4.3 The Rest of the Windmill Software Suite

The Windmill suite of software that comes with the 751 comprises: ConflML; SetupIML; the display and control panels—AnalogOut, DigitalOut & DDE Panel; Logger and Chart.

The display and control panels let you send data to, or display data from, any number of analogue and digital channels. Logger logs data to disk from up to 100 channels whilst Chart displays moving charts of data from up to 8 channels. For more channels, or different logging and charting speeds, simply run more instances of Logger and Chart. Full details of all these programs are given in their on-line Help files and the *Windmill User Manual*.

Should you need more sophisticated analysis or presentation, you can share data with other Windows applications using dynamic data exchange (DDE). For example, you can process data as it's collected using Microsoft Excel. See the DDE Panel Help file for more details.

There are many other optional programs in the Windmill range. **Graphics** lets you design and create your own Windmill displays—process mimics, wiring diagrams, bar charts, annunciator panels...whatever you wish. **Test-Seq** interprets a file of commands, and controls a test-rig accordingly. **Replay** replays a data file graphically. For details of other optional programs see <http://www.windmill.co.uk/programs.html>.

To be informed when new Windmill products are launched, be given data acquisition hints and tips, and read articles on measurement and control; subscribe to the free *Monitor* newsletter (ISSN 1472-0221) at <http://www.windmill.co.uk/newsletter.html>.

Index

- 590 screw terminal unit 2.16
- 590-8A current unit 2.16, 2.18
- 593 thermocouple unit 2.6, 2.18
- 594 bridge unit 2.7, 2.19, 2.21
- ac input 2.16
- analogue input 2.4
- auxiliary input 2.4, 2.6, 2.18
- biasing to ground 2.16
- bridge input 2.7, 2.19
- changing hardware settings 3.4
- cold junction 2.6, 2.18
- ConflML 2.4, 3.3-7
- contact closure 2.9, 2.16
- counter 2.11-12
- current input 2.18
- differential inputs 2.4
- digital input and output 2.9-10, 2.12, 3.2
- directory 3.6
- earth 2.4-5
- event counter 2.11
- filter 2.17-18
- high voltage 2.16
- id code 2.2, 3.4
- input 2.4, 2.9
- input filter 2.17-18
- input protection 2.16-18
- introduction 2.1
- isothermal box 2.18
- low pass filter 2.17
- new hardware 3.3
- noise 2.16-18
- output 2.9, 2.12
- plug 2.1
- power 2.1
- pressure transducer 2.19
- programmable resistance 2.16
- range 2.5
- resistance 2.16
- screw terminals 2.16
- SetupIML 4.2-6
- software
 - see Windmill software
- specifications 2.14-15
- strain gauge bridge 2.1, 2.7, 2.19
- temperature measurement 2.6, 2.18
- thermocouple 2.1, 2.6, 2.18
- totalise counter 2.11
- Windmill software 2.1, 2.4, 2.6, 2.18-20, 3.1-7
 - ConflML 3.3-7
 - earlier versions 3.2
 - installing 3.1-2
 - other modules 4.9
 - SetupIML 4.2-6
 - several copies 3.2
 - working directory 3.6